

From Blue-Ray, High Definition DVD to Near-Infrared Rays

An Advanced Optical Power Meter to Meet Next Generation Needs



A Truly Flexible Instrument Catering to Applications in R&D,

Optical Power Meter 3664 serves as a convenient tool for the R&D, manufacture and maintenance of equipment that incorporate laser light sources, such as DVD recorders, CD drives, copiers, and laser printers. With the addition of the Optical Sensors 9743 and 9743-10 specifically designed for blue-violet optical lasers that have low dependency on the incidence angle and flat wave sensitivity characteristics over its 405 nm bandwidth, the 3664 also proves to be a powerful instrument for inspecting next generation optical pickup drive devices.

Four Sensors to Accommodate All Types of Testing Needs


Choose from 4 different optical sensors according to wavelength and sensor structure to aptly meet your application requirements.

■ Two Different Bandwidths

9743/9743-10: optional sensor for blue-violet optical lasers only

Incorporated with characteristics that are ideal for measuring blue-violet lasers found in high-definition DVDs, Models 9743 and 9743-10 offer the following features:

- Planar wavelength sensitivity characteristics at the 405 nm bandwidth
- Low Incident Angle Dependency and Low Reflection

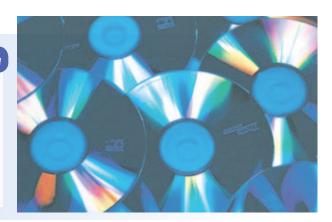



Test DVD and CD drives, Copiers, Printers and much, much more.

• Wide 320 to 1100 nm wavelength



■ Two Different Styles







Production and Maintenance of Laser- and LED-Dependent **Equipment**

Powerful Features

Superior cost performance

Guaranteed accuracy of ±5%, ideal for the production and testing of optical pickup devices (in combination with the Optical Sensor 9742/9742-10/9743/9743-10) Scaling functionality and USB compatibility all for one reasonable price.

Scaling

Adjust for sensitivity at the wavelength level, and easily integrate the 3664 into an inspection standard device.

■ Wavelength setting resolution up to 1 nm

Up to 10 wavelength memory presets can be configured, including the defaults for each optical sensor. The default wavelengths for each optional sensor are already stored in memory. Add up to 6 more custom wavelengths and call them up simply by pressing the λ key.

■ MAX/MIN/AVG measurement

Display maximum, minimum, and averages, as well as make relative measurements.



■ 2 Power settings

Compatible to both DC power and AC power. Choose a power method to suit your application.

Relative measurement

Display measurements as relative values, i.e., the difference from a set reference value. Load reference values from a measured value or define according to your requirements.

■ Upload data through the **USB** interface

Program your PC to download captured data, configure and even control the 3664 through the USB interface.

(USB driver software included)

System requirements

- OS: Windows 8/7/ Vista (32 bit/ 64 bit) Hardware requirements, such as those involving the CPU, RAM, and display monitor, shall conform to the system requirements of the OS.
- HDD space: 10 MB or more of free disk space
- Interface: USB Ver 1.1 or higher (connectable to only one PC)

Specifications

■ OPTICAL POWER METER 3664 Specifications

| Optical power measurement | | Units W/dBm |
|--|-----------------------------|--|
| | Range | Auto (manual settings available) |
| A | ccuracy | \pm 0.7 % (\pm 5 % when used with optional light sensor) |
| Ca | alibration | Resolution of 1 nm, automatic calibration of sensor, up to 10 |
| wa | avelength | wavelength presets available (including defaults for each sensor) |
| | Scaling | Configurable for each wavelength |
| | | Displays a measured value compared with a reference value |
| | otical loss | (Displayed value = measurement - reference) |
| mea | easurement | * Reference value can be based on a measurement, or input manually * Settings range: 0.001 nW to 1.9999 W (-90.00 dBm to 33.00 dBm) |
| | Display | 4 $\frac{1}{2}$ digits, up to 19999. Display resolution: 0.01 dBm/0.01 dB |
| | | Units: nW/µW/mW/dBm/dB |
| Measurement display Wavelength display | | 4 digits, unit: nm |
| Display refresh rate | | Approx. 330 ms |
| | , | 11 |
| | /MIN display | Displays MAX/MIN during measurement |
| AV | E display | Moving average, average count configurable from 2 to 100 |
| A | nalog Output | According to optional Optical Sensor in use |
| Ana | | 9742/9742-10: 1V at sensor correction input 9743/9743-10: 0.7V at sensor correction input |
| Outn | ut resistance | 50Ω |
| | | |
| Outp | tput connector Interface | φ3.5 mini jack |
| Ir | | USB Ver1.1 Output of measurement data, configuration and control supported |
| Inclu | ded features | Auto power save, configuration backup, battery check |
| | | Safety standard: EN61010 |
| | oplicable andards | EMC:EN61326, EM61000-3-2, EM61000-3-3 |
| | Power | LR6 (AA) Alkaline battery ×4, AC adapter (9445-02/9445-03) |
| | | 1.6 VA |
| Max. rated power | | |
| Ope | perating time | 60 hours (when using battery, continuous use with 9742 optical sensor as correction input) |
| Operating conditions | | 0°C to -40°C (32°F to 104°F), up to 80 % rh (no condensation) |
| Stora | ige conditions | -10°C to 50°C (14°F to 122°F), up to 80 % rh (no condensation) |
| Operating environment | | Indoor, up to elevation of 2000 meters (6562 ft) |
| | noiono and mass | 85W × 160H × 35D mm (3.35W × 6.30H × 1.38 D inches) |
| Dimensions and mass | | (excluding protrusions), 270g (9.5 oz.) (without batteries) |

Accuracy: 23°±5°C (73°±9°F), less than 80% Accuracy guaranteed for 1 year, Post-adjustment accuracy guaranteed for 1 y

OPTICAL SENSOR Specifications

0749/0749 10

| 9742/ 9742-10 | | |
|-----------------------------------|--|--|
| Measured wavelength | 320 to 1100 nm | |
| Measured power | -59 dBm to +17 dBm (correction wavelength) | |
| Maximum rated value | 50 mW (+17 dBm) (under direct lighting) | |
| Receiving element | Si Photo diode | |
| Receptor size | Approx. 9.6 mm × 9.6 mm (0.38" × 0.38") | |
| Accuracy | \pm 4.3 % (± 5 % when used with Optical power meter 3664) [Correction conditions] correction wavelength 633 nm, correction power 100 μ W, when ø approx. 2 mm (0.08") parallel beam strikes perpendicular to center of optical sensor, CW light | |
| Wavelength configuration defaults | 633 nm, 635 nm, 650 nm, 780 nm | |
| Operating conditions | | |
| Storage conditions | -10 °C to 50 °C (14°F to 122°F), up to 80 % rh (no condensation) | |
| Operating environment | Indoor, elevation up to 2000 meters (6562 ft) | |
| Mass | 100 g (3.5 oz.) | |
| 9743/ 9743-10 | | |
| Measured wavelength | | |
| Measured power | -50 dBm to + 20 dBm (correction wavelength) | |
| Maximum rated value | 100 mW (+20 dBm) (under direct lighting) | |
| Receiving element | | |
| Receptor size | 10 mm × 10 mm (0.39" × 0.39") | |
| Accuracy | $\pm 4.3\%$ (±5% when used with Optical power meter 3664) [Correction conditions] correction wavelength 405 nm (Using a 405 ± 5 nm wavelength as the reference wavelength), correction power 100 μ W, when ϕ 1.5 mm (0.06") parallel beam strikes perpendicular to center of optical sensor | |
| Wavelength configuration defaults | 400 nm, 403 nm, 405 nm, 408 nm | |
| Operating conditions | 0 °C to -40 °C (32°F to 104°F), up to 80 % rh (no condensation) | |
| Storage conditions | -10 °C to 50 °C (14°F to 122°F), up to 80 % rh (no condensation) | |
| Operating environment | Indoor, elevation up to 2000 meters (6562 ft) | |
| Mass | [9743]: 110 g (3.9 oz.) [9743-10]: 100 g (3.5 oz.) | |

Please refer to the external diagrams depicted on the next page for the exact dimensions of the optical sensors.

OPTICAL SENSOR exterior view

■ OPTICAL SENSOR 9742 ■ OPTICAL SENSOR 9742-10 **■ OPTICAL SENSOR 9743** ■ OPTICAL SENSOR 9743-10 φ2.7×2 Ŧ 160 160 Cable Cable 180 180 72.5 84 84 Lagged 40 18 40 30 (can be secured with screws) (can be secured with screws) (can be secured with screws)

Model: OPTICAL POWER METER 3664


Model No. (Order Code) (Note)

3664

Accessories: AC adapter 9445-02 (for USA) ×1 or 9445-03 (for EU) ×1, Output cord L9094 ×1, Driver software (CD-R) ×1, LR6 (AA) alkaline batteries ×4, USB cable ×1, Strap ×1, Instruction manual ×1

The OPTICAL POWER METER 3664 must be used in conjunction with either the Optical sensor 9742 or 9742-10 or 9743 or 9743-10, sold separately.



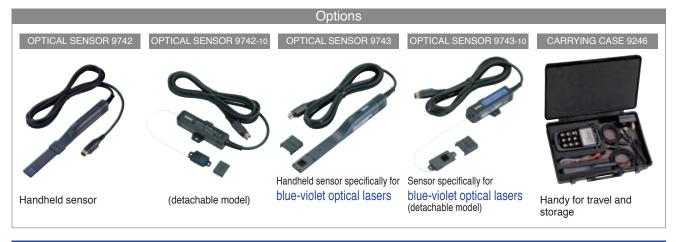

Optical Power Meters installed with firmware version 1.01 or earlier must be updated to support compatibility with the new Optical Sensor 9743/ 9743-10

When using Model 9743/9743-10 with an earlier version of Model 3664, upon connecting the sensor to the optical power meter and turning on the power, "Err1" will be displayed.

To update your Model 3664, please visit our website at www.hioki.com to download the latest firmware, "HIOKI 3664 Setup Utility" and refer to the instructions provided to learn how to set up the proprietary data for the new optical sensor.



HEADQUARTERS

81 Koizumi, Ueda, Nagano, 386-1192, Japan TEL +81-268-28-0562 FAX +81-268-28-0568 http://www.hioki.com/E-mail: os-com@hioki.co.jp

HIOKI (Shanghai) SALES & TRADING CO., LTD. TEL +86-21-63910090 FAX +86-21-63910360

http://www.hioki.cn / E-mail: info@hioki.com.cn

HIOKI SINGAPORE PTE. LTD. TEL +65-6634-7677 FAX +65-6634-7477 E-mail: info-sg@hioki.com.sg

HIOKI KOREA CO., LTD.

TEL +82-2-2183-8847 FAX +82-2-2183-3360 E-mail: info-kr@hioki.co.jp

HIOKI USA CORPORATION

TEL +1-609-409-9109 FAX +1-609-409-9108 http://www.hiokiusa.com / E-mail: hioki@hiokiusa.com

Note: Company names and Product names appearing in this catalog are trademarks or registered trademarks of various companies.

DISTRIBUTED BY